

Tissue Culture Facility Protocol

- No bacteria culture should be brought to the facility.
- The incubator in the facility should be used as temporary storage only. Clean up any spill in the incubator with ethanol. If media spilled into the water tray, clean up the tray and fill with clean dI water.
- Disinfect all liquid waste (including the vacuum flask) with bleach and pour down the drain.
- Dispose all bio-waste in the red trash. Anything left in the centrifuge or on the bench top will be tossed as waste.
- When you are done
 - Wipe off the biosafety cabinet with ethanol,
 - Turn off the light, vacuum and receptacle,
 - Leave the air flow on and the cabinet shield open.
 - Turn off the water bath, centrifuge and the microscope.
 - Make sure the door of the CO₂ incubator is closed all the way.
 - Wash your hands.
 - Turn off the light.