

Announcements

- Mid-term feedback and compoment grades will be up on website by next week
- Chapter 9 Laboratory Report Template will be posted by Wednesday (3/26)
- Chapter 9 Raw Data file and Laboratory Report due April 3 – 8 (>1 week; ~2 weeks from when you finish the lab)

Announce

Concepts

Procedure

Hazards

Tips

Clarification

End

Chapter 9 Calculations: Comparing two methods of protein labeling

- Figure 1: BCA standard curve
- * • Table 1: Protein Concentration (μM) for both **DTNB** and **F5M** reactions
- Figure 2: Effect of **DTNB** on KHK rates.
- * • Table 2: Effect of **DTNB** on KHK specific activity
- Figure 3: Effect on Coupled assay of [PK].
- Figure 4: Plot of all **DTNB** reactions.
- * • Table 3: Calculate μM Cys from **DTNB** data.
- Figure 5: **F5M** standard curve
- * • Table 4: Calculate μM Cys from **F5M** data
- * • Table 5: Calculate the number of Cys per KHK molecule by these two different labeling methods
- Table 6: Show calculation of error propagation

Announce

Concepts

Procedure

Hazards

Tips

Clarification

End

Error Propagation in Chapter 9

Purpose:

- To account for uncertainty in measured data which will carry over into subsequent steps and assays
- Often denoted as $X \pm x$ where,
 - X = data measurements
 - x = error values
 - $(x/X)100$ = (fractional error) percent error

You will need to take error propagation into account for ALL steps in:

Chapter 9A

- Measuring the effect of DTNB on KHK activity (Tbl2)
- Measuring the amount of DTNB that reacted with Cys residues in KHK (Tbl3)

Chapter 9BE

- Measuring RFUs vs. μM (or moles) of F5M in standard curve (Fig5)
- Measuring μM (or moles) of F5M labeled on KHK (Tbl4)
- Measuring μM (or moles) of KHK in DTNB & F5M reactions from BCA assay (Tbl1)
- Measuring A_{562} vs. μM (or moles) BSA in BCA assay in standard curve (Fig1)
- Calculate number of Cys by two methods (Tbl 5&6)

Announce Concepts Procedure Hazards Tips Clarification End

Overview of Chapter 9 Experiment

Determination of number of Cys residues in KHK

Table 3

Table 4

Table 5

DTNB:

$$\frac{A_{412} \times 0.00025 \text{ L}}{("14,000" \text{ M}^{-1}\text{cm}^{-1})(\text{"x cm"})} = \text{number of moles Cys in KHK in the 250 } \mu\text{L reaction (D and F)}$$

$$\frac{\text{mg of KHK protein in the 250 } \mu\text{L reaction} \times 0.001 \text{ g/mg}}{\text{MW of KHK subunit "34,000" g/mole}} = \frac{\text{moles Cys}}{\text{moles KHK}}$$

F5M:

Concentration of F5M labeled to KHK $\times \sim 0.000035 \text{ L}$ = number of moles Cys in KHK in the 35 μL reaction (A and B)

$$\frac{\text{mg of KHK protein in the eluted } \sim 35 \mu\text{L reaction} \times 0.001 \text{ g/mg}}{\text{MW of KHK subunit "34,000" g/mole}} = \frac{\text{moles Cys}}{\text{moles KHK}}$$

Need to account for the various dilutions you subjected your samples to, e.g., 10x dilution

Can be simplified if use μM as the reaction volumes cancel out

- Compare the two numbers and determine if they agree with each other or not
- Are the values statistically different? – *Important to do error propagation!*

Table 6

Announce Concepts Procedure Hazards Tips Clarification End

Error Propagation formulas

- Follow these formulas for propagating data measurements with error

Addition: $(X \pm x) + (Y \pm y) = (X + Y) \pm \sqrt{x^2 + y^2}$

Subtraction: $(X \pm x) - (Y \pm y) = (X - Y) \pm \sqrt{x^2 + y^2}$

Multiplication: $(X \pm x) \times (Y \pm y) = (X \times Y) \pm (X \times Y) \sqrt{\left(\frac{x}{X}\right)^2 + \left(\frac{y}{Y}\right)^2}$

Division: $(X \pm x) \div (Y \pm y) = (X \div Y) \pm (X \div Y) \sqrt{\left(\frac{x}{X}\right)^2 + \left(\frac{y}{Y}\right)^2}$

Calculated result \swarrow \nwarrow Calculated error

Announce

Concepts

Procedure

Hazards

Tips

Clarification

End

Make your own error propagation calculator using excel!

Error Propagation Calculator																							
Values below are entered as: (A \pm a) <function> (B \pm b) = (C \pm c)																							
A	\pm	a																					
10.00000	\pm	0.01000																					
B	\pm	b																					
5.00000	\pm	0.05000																					
<table border="1"> <thead> <tr> <th></th><th>C</th><th>\pm</th><th>c</th></tr> </thead> <tbody> <tr> <td>Addition</td><td>15.00000</td><td>\pm</td><td>0.0509902</td></tr> <tr> <td>Subtraction</td><td>5.00000</td><td>\pm</td><td>0.0509902</td></tr> <tr> <td>Multiplication</td><td>50.00000</td><td>\pm</td><td>0.5024938</td></tr> <tr> <td>Division</td><td>2.00000</td><td>\pm</td><td>0.0200998</td></tr> </tbody> </table>					C	\pm	c	Addition	15.00000	\pm	0.0509902	Subtraction	5.00000	\pm	0.0509902	Multiplication	50.00000	\pm	0.5024938	Division	2.00000	\pm	0.0200998
	C	\pm	c																				
Addition	15.00000	\pm	0.0509902																				
Subtraction	5.00000	\pm	0.0509902																				
Multiplication	50.00000	\pm	0.5024938																				
Division	2.00000	\pm	0.0200998																				

Announce

Concepts

Procedure

Hazards

Tips

Clarification


End

Error Propagation from standard curves that are not absolutely linear: BCA & F5M

Fig. 1
Table 1
Fig. 5
Table 4

Method 1:

- Repeat assay on several identical aliquots of your proteins. Use average and standard deviation of measured values from standard curve. *(This is what you will be doing in 9C!)*

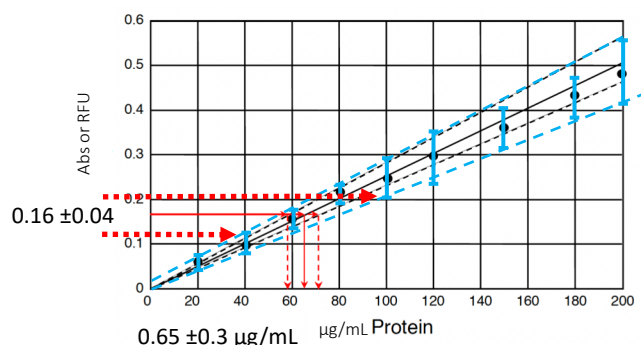

Method 2:

- ❖ Use LINEST function in Excel based on your standard curve

<https://www.youtube.com/watch?v=BvQ11wtKeCg>

Announce Concepts Procedure Hazards Tips Clarification End

Error Propagation from BCA assay to determine protein concentration

To calculate moles of KHK in your DTNB and F5M reactions:

- Find KHK concentration of DTNB assay / F5M labeling reaction by making use of the BSA standard curve you created and measured triplicate absorbance values. Take the average and standard deviation in your raw data workbook.* Use to get concentration in the 25 µL of sample you measured in µg/mL.
- Since we have the same volumes from the assay as in the standard curve, and they are directly from the reaction without dilution, we can convert µg/mL to µM using the molecular weight of KHK (**34,000 g/mol**; the error in MW is so low (<0.01%), it can be ignored)

➤ Example: $[(275 \pm 25 \text{ µg/mL})(10^{-6} \text{ g/µg})(10^3 \text{ mL/L})] \div [34000 \text{ g/mol}]$
 $= 8.1 \pm 0.7 \times 10^{-6} \text{ mol/L} \times 10^6 \text{ µmol/mol} = 8.1 \pm 0.7 \text{ µM}$

➤ [The easy calculation is $\text{µg/mL} \div 34 = \text{µM}$]

* STDEV uses the following formula:

$$\sqrt{\frac{\sum (x - \bar{x})^2}{(n-1)}}$$

where \bar{x} is the sample mean AVERAGE(number1,number2,...) and n is the sample size.

Announce Concepts Procedure Hazards Tips Clarification End

How to convert DTNB signal to moles of Cysteine residues?

Fig. 1
Table 1
Fig. 5
Table 4

To calculate μM –SH groups from DTNB modification:

- Calculate the average **net** absorbance by subtracting the background
 - Propagate the error by using subtraction equation
- Use **Beer's Law** $\rightarrow A = \epsilon lc$
 - End-point average net absorbance = extinction coefficient x path length x concentration
 - Extinction coefficient @ 412nm for DTNB = $14,000 \text{ M}^{-1}\text{cm}^{-1}$ (or $0.014 \mu\text{M}^{-1}\text{cm}^{-1}$)
 - Path length of 250 μL sample in 96-well plate is 1 cm
 - Solve for concentration **in μM** by taking the highest absorbance reading of your sample after 1 h run time, or the last from your plot
- Calculate Error in your determination
 - Propagating the error by using the same fractional error (since ϵ & l have negligible errors, OR
 - Repeat entire calculation of 3 trials and calculate the standard deviation

Subtraction: $(X \pm x) - (Y \pm y) = (X - Y) \pm \sqrt{x^2 + y^2}$

Announce

Concepts

Procedure

Hazards

Tips

Clarification

End

How to convert F5M RFU signal to moles of Cysteine residues?

Fig. 5
Table 4

To calculate μM –SH groups from F5M reactions:

- Use **standard curve** which relates concentration of F5M with RFU
 - Calculate **net** RFU and error for each sample
 - Propagate the error by using subtraction equation
- Use graph to find concentration and error in μM from your sample **net** average RFU values and standard deviations

Subtraction: $(X \pm x) - (Y \pm y) = (X - Y) \pm \sqrt{x^2 + y^2}$

Chapter 10

LIPIDS AND MEMBRANES

Chapter 10: Overview

- **Week 1 – 10A:**
 - Isolate and wash erythrocytes and plasma membrane fractions (washing & centrifugation)
- **Week 2 – 10BC:**
 - Determine yield of membrane, lipids, and composition of lipids (TLC & cholesterol assay)
- **Week 3 – 10D:**
 - Separation of proteins by SDS-PAGE and immunoblotting
- **Week 4 – 10E:**
 - Visualization of SDS-PAGE and immunoblot results (i.e. Western Blot analysis)

Chapter 10A: Lipids and Membranes

Objectives

- Understand lipid membrane composition, structure, and function
- To isolate red blood cells from bovine (*Bos taurus*) blood
- Prepare washed and unwashed membrane fractions

Procedures

- **Isolate erythrocytes** from fresh bovine blood by **centrifugation**
- **Lyse erythrocytes** via **osmotic shock using low-salt buffer**
- To fractionate membrane preparations containing peripheral bound proteins away from integral membrane proteins
- **Lyophilize** (freeze dry) membrane preparations for next week's lipid extraction

Announce

Concepts

Procedure

Hazards


Tips

Clarification

End

Plasma Membrane

- Physical barrier between cell and external environment
- Separates organelles and subcellular structures
- Composed of lipids, proteins, and carbohydrates
- Exists in a **fluid-mosaic model**
 - Lipids expose **hydrophilic** head to aqueous environment
 - **Hydrophobic** lipid tails are internal to membrane
 - Lipid-bilayer embedded with proteins

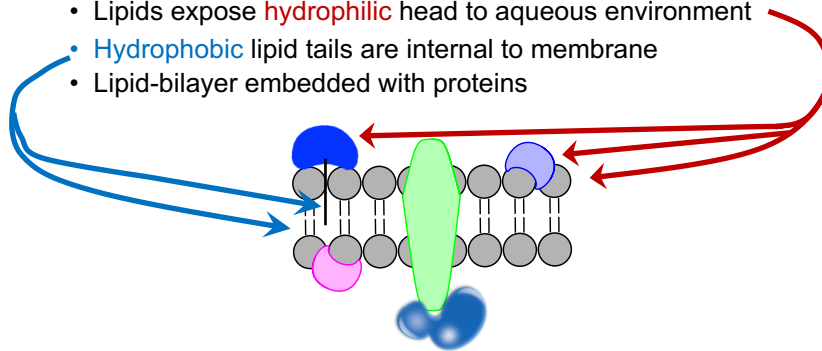

Announce

Concepts

Procedure


Hazards

Tips

Clarification

End

Plasma Membrane

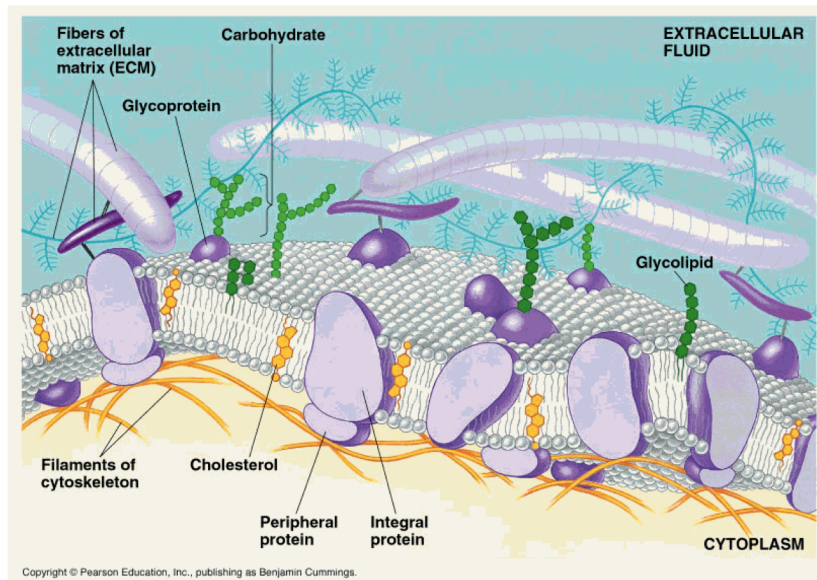

Announce

Concepts

Procedure

Hazards

Tips

Clarification

End

Erythrocyte Membranes

- **Red Blood Cells – Erythrocytes**

- Human Erythrocyte Membrane:

- 50% protein
- 40% lipid
- 10% carbohydrate



- Isolated as “ghosts” –remains after removing hemoglobin and intracellular components
- Isolated from whole blood extract through **cell lysis and centrifugation**
- Do not contain other organelles – easier sample preparation

Announce

Concepts

Procedure

Hazards

Tips

Clarification

End

Isolation of Membranes

- Red Blood Cells **ISOLATED** from blood with low-speed centrifugation using Phosphate Buffered Saline (**PBS**)
 - Platelets and lymphocytes (white blood cells) float to top, RBCs to bottom
- RBCs are **LYSED** using Hypotonic Phosphate Buffered Saline (**HPBS**)
 - Low osmotic pressure in HPBS buffer – cells swell and rupture due to hypotonic solution
 - Only plasma membrane remains, because no organelles in RBCs

Announce

Concepts

Procedure

Hazards

Tips


Clarification

End

How does HPBS rupture erythrocytes?

Osmolysis

If **EXTRACELLULAR ENVIRONMENT** is:

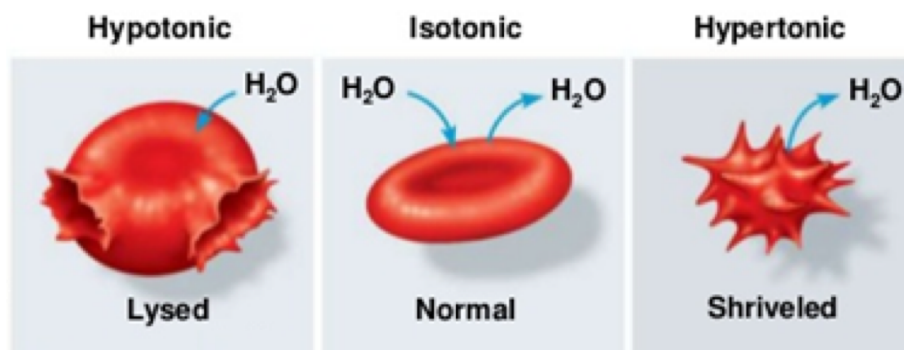

Announce

Concepts

Procedure

Hazards

Tips

Clarification

End

Isolation of Membrane Components

- Membrane components can be dissolved and separated by various solvents:
 - **Lipids** – organic solvent mixtures
 - **Carbohydrates** – typically found associated with proteins
 - **Proteins**
 - **Integral** – treating with detergents
 - **Peripheral** – washing with a high-salt buffer

Announce

Concepts

Procedure

Hazards


Tips

Clarification

End

Isolation of Membrane Components

- **3 Fractions:**
- Unwashed Membrane
 - **Peripheral and Integral**
- Washed Membrane
 - **Integral**
- Washed Supernatant Fraction
 - **Peripheral**

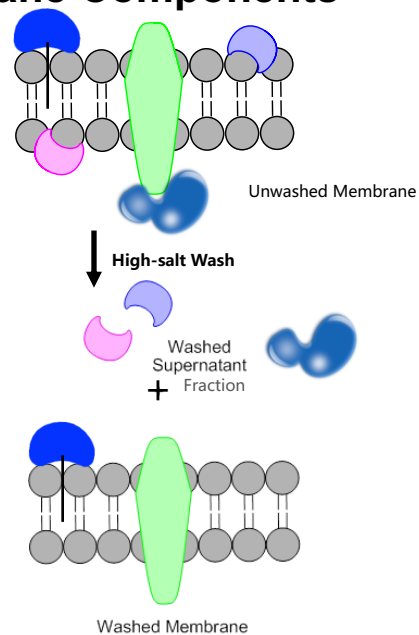

Announce

Concepts

Procedure

Hazards

Tips

Clarification


End

Isolation of Membrane Components

- **3 Fractions:**
 - Unwashed Membrane
 - **Peripheral and Integral**
 - Washed Membrane
 - **Integral**
 - Washed Supernatant Fraction
 - **Peripheral**
- **Lyophilization**
 - Freeze drying of membrane preparations
 - Removes water by sublimation
 - Leaves all solids
- **Solids will be used to extract lipids with organic solvents**



LYOPHILIZATION

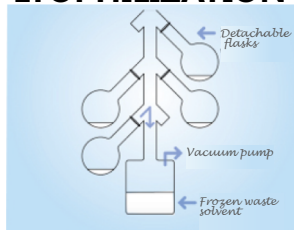

Freeze dried raspberries



Announce

Concepts

Procedure

Hazards



Tips

Clarification

End

Part I: Isolating erythrocytes

Keep everything on ice!

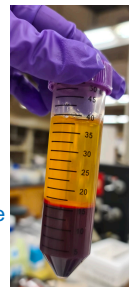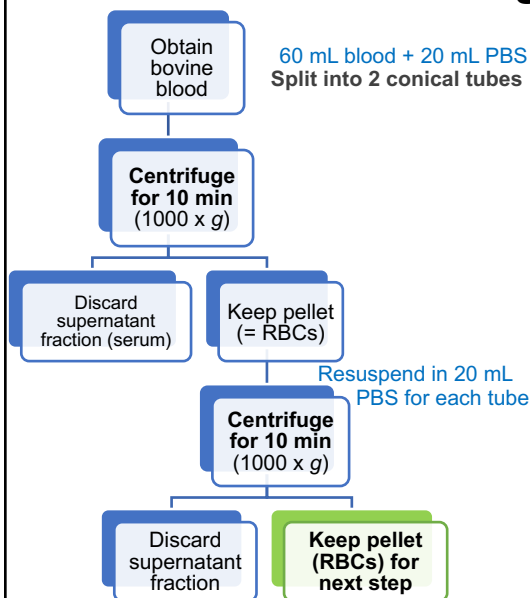

- Use swinging-bucket centrifuge rotors for 50 ml conical tubes
- **Make sure to balance tubes!**
- **Make sure to use ONLY PBS – NOT HPBS!!!**
(Remember why??)
- Some steps may be done in advance by TFs

Announce

Concepts

Procedure


Hazards

Tips


Clarification

End

Part II: Isolating the plasma membrane

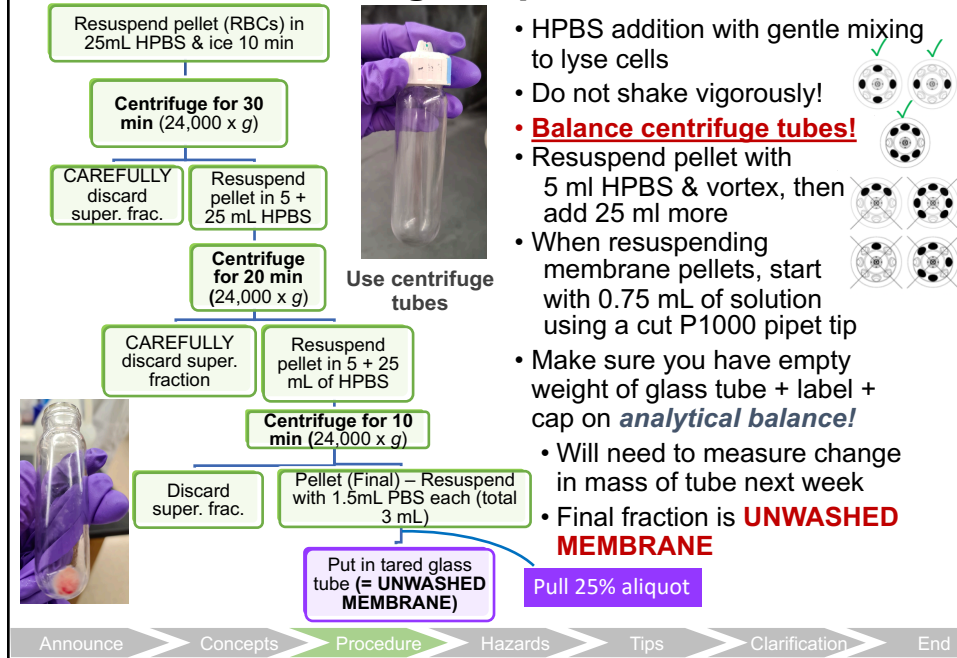

Part III: Washing the membrane

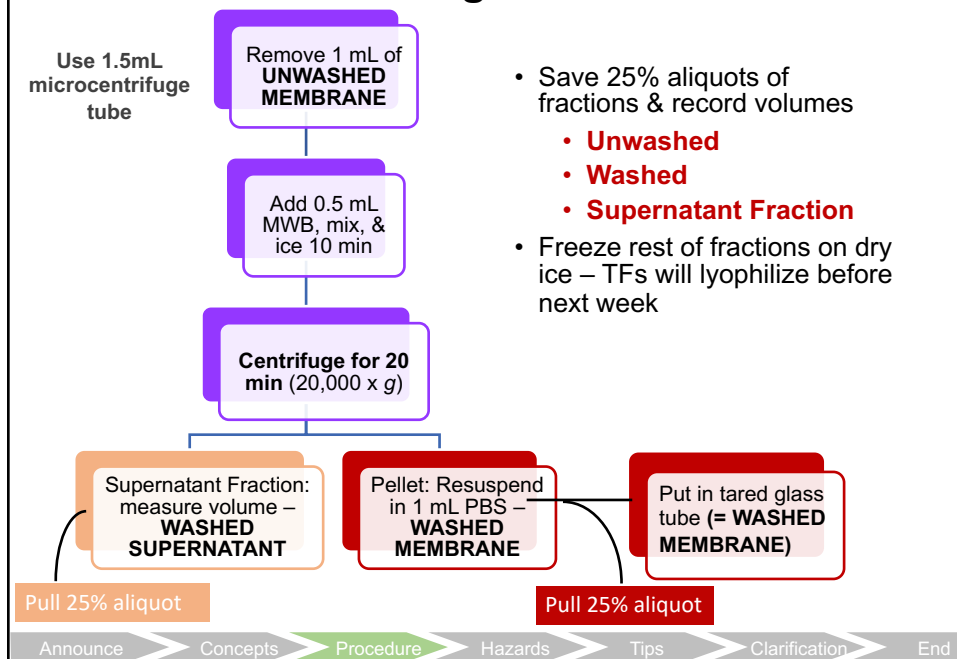

CH 10A waste disposal

- Blood waste can go into a **separated** nearby collection carboy in the fume hood
- All other reagents are safe to dispose into the main steam waste carboy

Announce > Concepts > Procedure > **Hazards** > Tips > Clarification > End

Chapter 10a Lab Tips

Work efficiently between centrifugation steps! (there are 1 h 40 min of spin time)

Reminders for accurate taring

1. Pre-tape and write **ALL** information on tubes **and** caps **before** taring
2. Stick to one analytical balance for the remainder of Ch 10
3. Do not shake or move benches when weighing out tubes
4. Wait 5-10 seconds after shutting glass screen to record mass

At the end of lab you will turn in the following:

In tared labeled glass tubes (+caps) --

1. 75% unwashed membrane
2. 75% washed membrane

In 1.5 mL eppendorf tubes --


3. 25% aliquot of unwashed membrane
4. 25% aliquot of washed membrane
5. 75% aliquot of washed supernatant fraction
6. 25% aliquot of washed supernatant fraction

Don't forget to ALWAYS record your measurements of mass and volume when directed to do so in procedures

Announce > Concepts > Procedure > Hazards > **Tips** > Clarification > End

Chapter 10A clarification

- First step involves preparing 2 tubes of 30 mL bovine blood + 10 mL PBS. Subsequent volume mentioned in the manual refers to each of the tubes

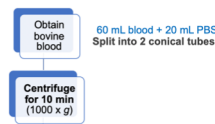

- Combine the two tubes of 1.5 mL each of the final pellet → UNWASHED membrane fraction



- For the step involving removal of 1.0 mL of unwashed membrane preparation to 1.5 mL microcentrifuge tube, use the P1000 pipets. It should fit.
- Remember to balance centrifuge tubes for ALL centrifugation steps

Announce > Concepts > Procedure > Hazards > Tips > **Clarification** > End


Chapter 10A

Before the lab period, you should have:

- ✓ Completed your Pre-lab Write-up and submit on Gradescope
 - ✓ Title, purpose and procedures
 - ✓ Remember to include:
 - ✓ Flowchart for your membrane isolation procedure
 - ✓ Note masses to record
 - ✓ Masses for **tared** Pyrex **tubes & caps** for 75% washed & 75% unwashed membrane
 - ✓ Note **volumes** for all six fractions
 - ✓ (25% & 75%) x (Unwashed membrane, washed membrane, and washed supernatant fraction)

At the end of lab, you should have:

- ✓ Turned in your **six** aliquots to your TFs
 - ✓ Unwashed and washed membrane glass tubes, "parafilm"
 - ✓ Four eppendorf tubes and glass tube caps in a conical tube
- ✓ Washed your high-speed centrifuge tubes → Clean → (do not throw them away in the trash)

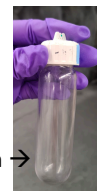

Announce > Concepts > Procedure > Hazards > Tips > **Clarification** > End

Questions?

Activity & Discussion Quiz