

Announcements

- Chapter 8 Week 1 by EOD Friday
- Chapter 8 Lab report template released Wed, Feb 11
- Chapter 7 Lab report grades will be returned by the Monday after the B5 section turns their reports in
- Lab feedback

Announcements

- Dr. Tolan here later to discuss Creativity Project

Chapter 8AB: In vitro transcription and translation

Objectives

- To use an in vitro transcription and translation system to synthesize proteins from genes cloned into a plasmid.
- Introduction to fluorescence

Procedures

- Use a **combined RNA polymerase transcription system** and a **wheat germ lysate-translation system** to ***express and synthesize proteins*** from cloned genes
- Use SDS-PAGE analysis to analyze protein expression

Transcription and Protein Synthesis



Week 1:

Add DNA template and perform transcription/translation using TNT® Wheat Germ Extract.



Week 2:

Separate translation products by SDS-PAGE



Fluorescent Imaging

Central dogma of molecular biology

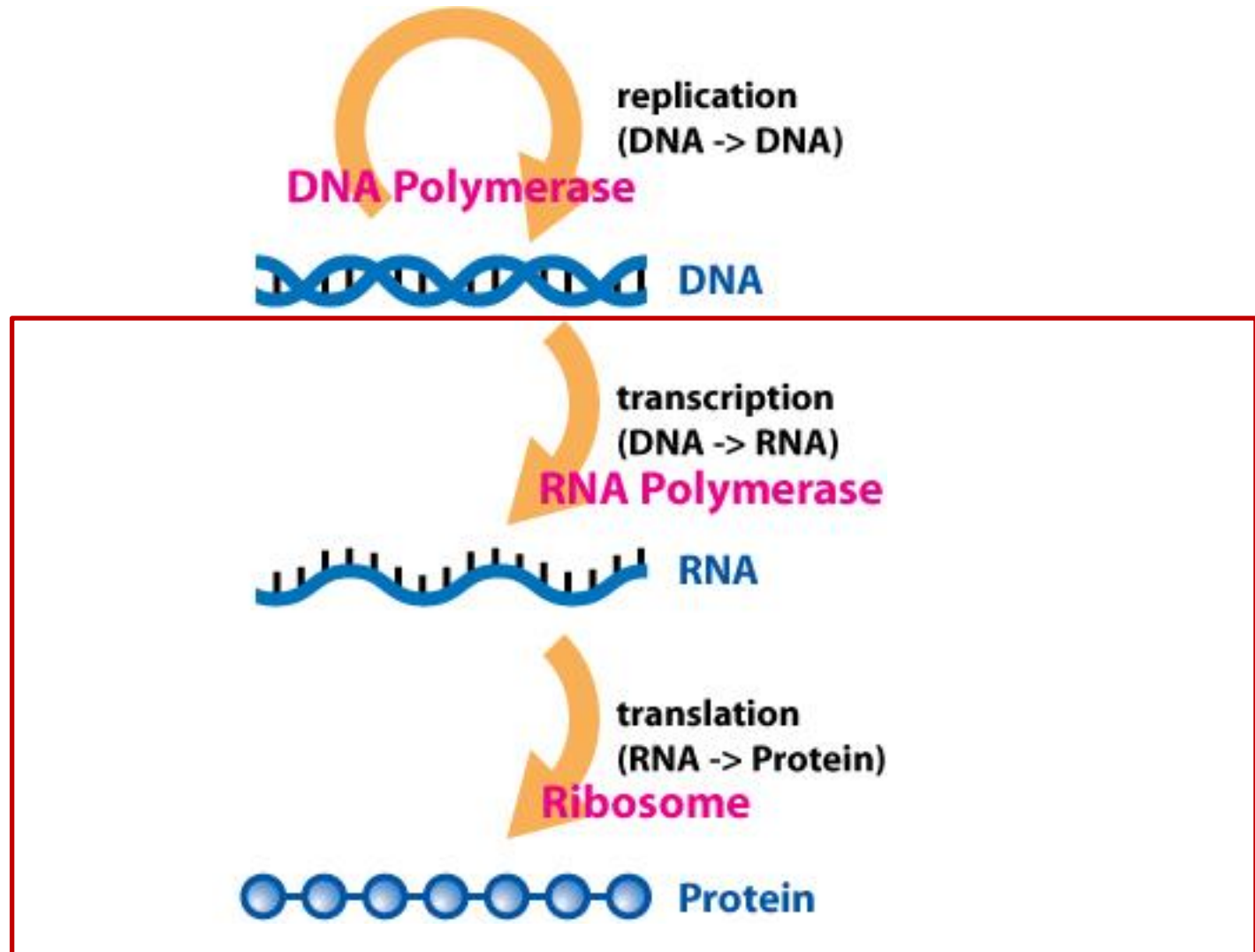


Photo: Wikipedia (2008).

In vitro transcription/translation

- **DNA systems:** coupled transcription/translation
- **RNA systems:** translation only

- Chapter 8 will use a **DNA system** (plasmid DNA > RNA > Protein)
 - Components in system include:
 - SP6 polymerase
 - NTPs (ATP, GTP, CTP, UTP)
 - Amino acids
 - Wheat Germ Extract
 - ❖ Ribosome and other translation factors
 - ❖ tRNAs
 - ❖ GTP (source of translation energy)
 - ❖ Fluortect Green_{Lys}
 - Students will supply the plasmid DNA


Recall from Chapter 6...

- Isolated pGEM3-Rel and pGEM4-Rel from *E. coli* using mini-prep
- Each plasmid contains:
 - **SP6 promoter**
 - **Target gene - Rel**
 - **Ampicillin resistance gene**
 - **Origin of replication**
 - **Restriction enzyme recognition sites**
- Transcription only starts at sites with specific promoter sequences
- Incorporated just upstream of cloning sites of any opening reading frame

Plasmid maps Figure 6-5, p. 209

Recall from Chapter 6...


- Isolated pGEM3-Rel and pGEM4-Rel from *E. coli* using mini-prep
- Each plasmid contains:
 - **SP6 promoter**
 - **Target gene - Rel**
 - **Ampicillin resistance gene**
 - **Origin of replication**
 - **Restriction enzyme recognition sites**
- Transcription only starts at sites with specific promoter sequences
- Incorporated just upstream of cloning sites of any opening reading frame

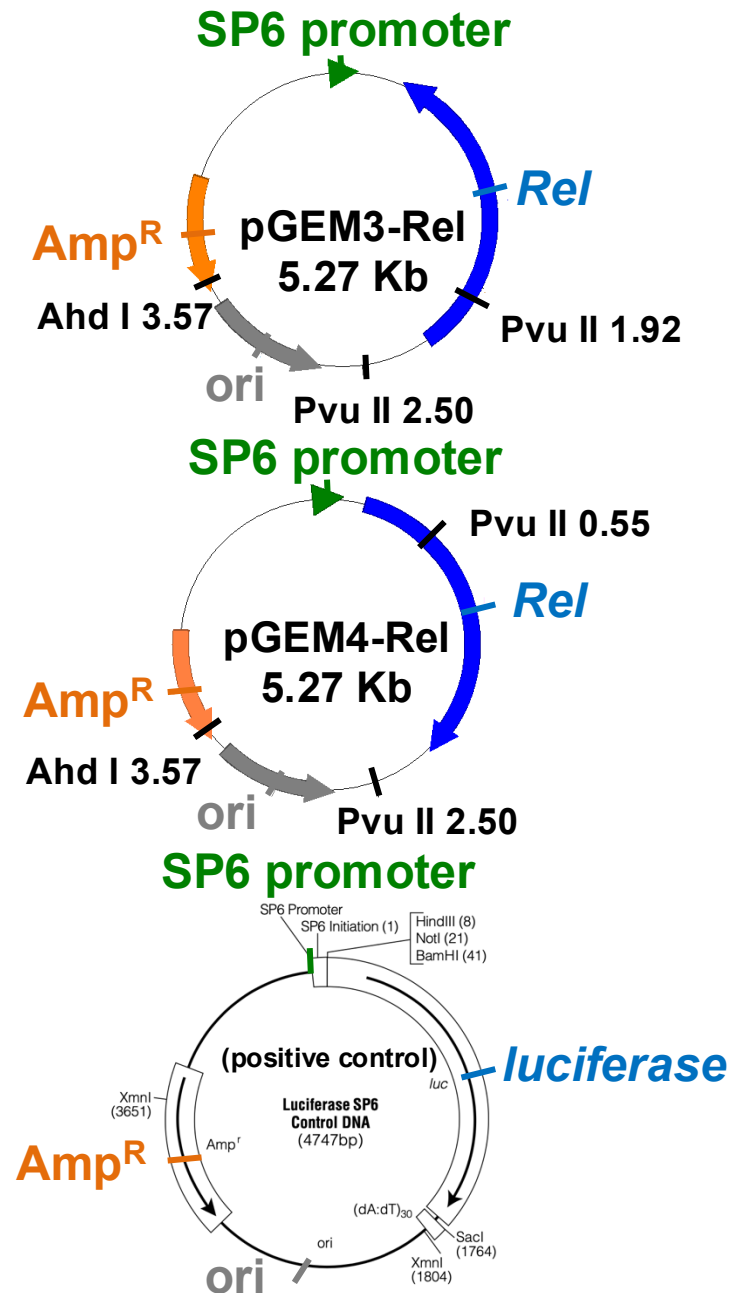

Plasmid maps Figure 6-5, p. 209

How will we know if it worked?

Use a
fluorescently
labeled
amino acid!

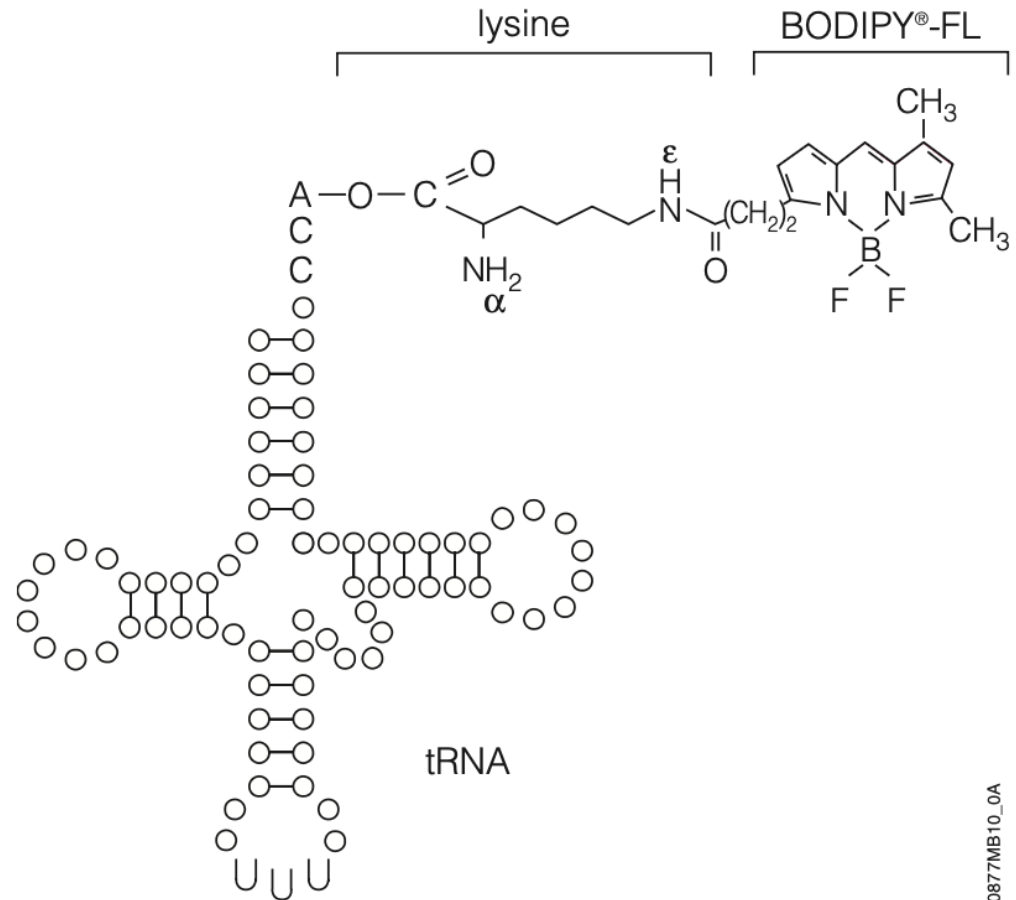

0877MB10_0A

Fluorescence spectroscopy

- Molecules that are capable of fluorescence are known as fluorophores
- Fluorophores absorb light energy at particular wavelengths and energy levels (absorption) and reach an excited state (excitation)






- It then emits light at longer wavelengths at a lower energy level (emission)



- Net reaction: $M + h\nu \rightarrow M^* \rightarrow M + h\nu'$

$$h\nu > h\nu'$$

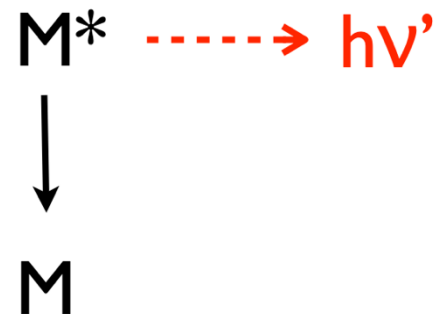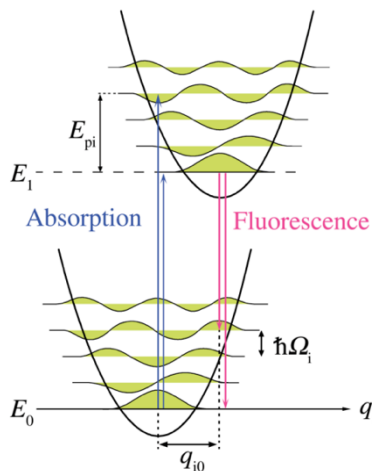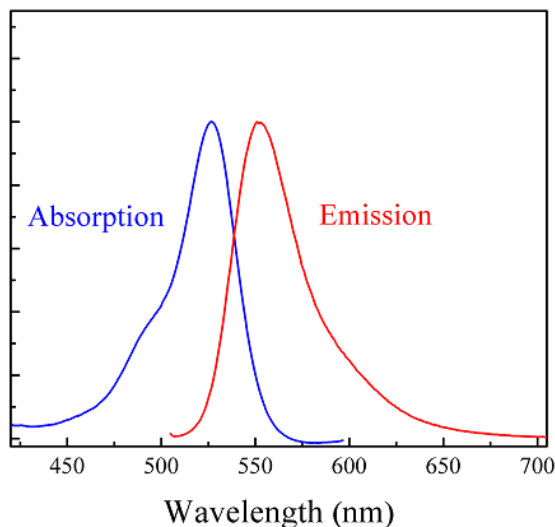

<https://www.chem.uci.edu/~unicorn/M3LC/handouts/Week2/SpectroscopyHandout.pdf>

Molecules and materials can be intrinsically fluorescent or synthesized

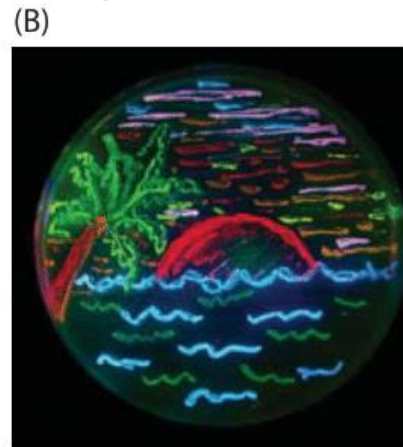
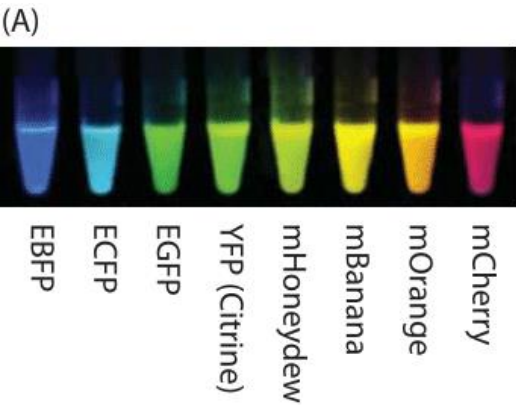


Illustration of some of the palette of fluorescent proteins that has revolutionized cell biology.




(A) Fluorescent proteins spanning a range of excitation and emission wavelengths. (B) Illustration of a petri dish with bacteria harboring eight different colors of fluorescent protein and used to “paint” an idyllic beach scene.

(Adapted from: R. Y. Tsien, Nobel lecture, *Integr. Biol.*, 2, 77-93, (2010).)



Examples of fluorescent molecules

- Amino acids (Trp, Phe, Tyr)
- Base pair derivatives (2-AP, 3-MI, 6-MI, 6-MAP, pyrrolo-C, tC)
- Chlorophylls
- Fluorescent Proteins (FPs)
- Organic dyes (fluorescein, rhodamine, N-aminocoumarins and derivatives of these)
- Rare earth elements (lanthanides)
- Quantum dots
- And many more...

Basic Flowchart of Methods

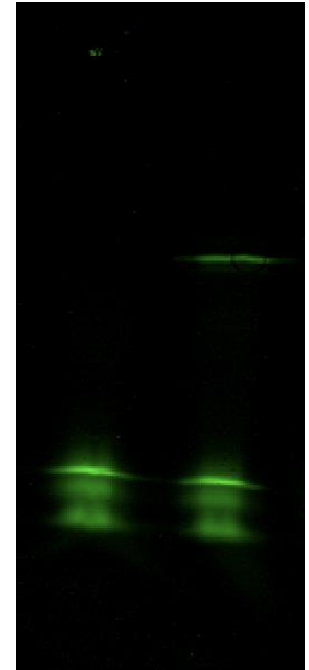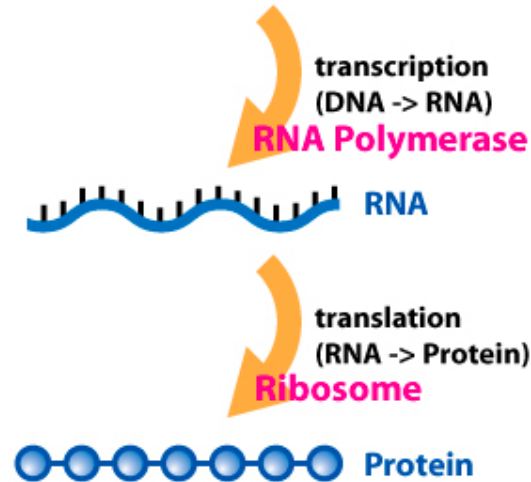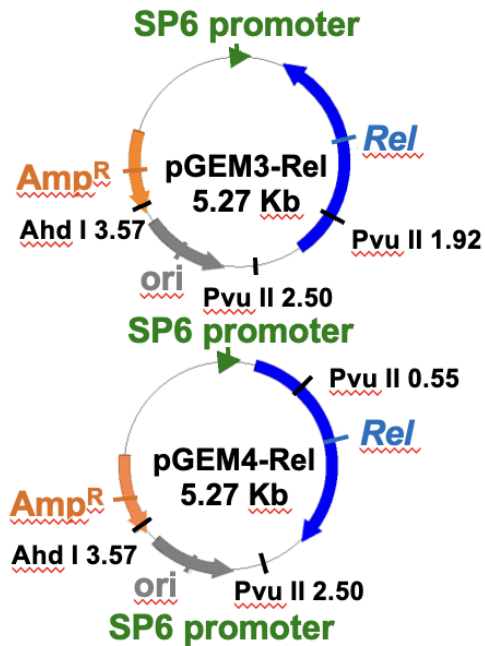

Intro to BODIPY-FL

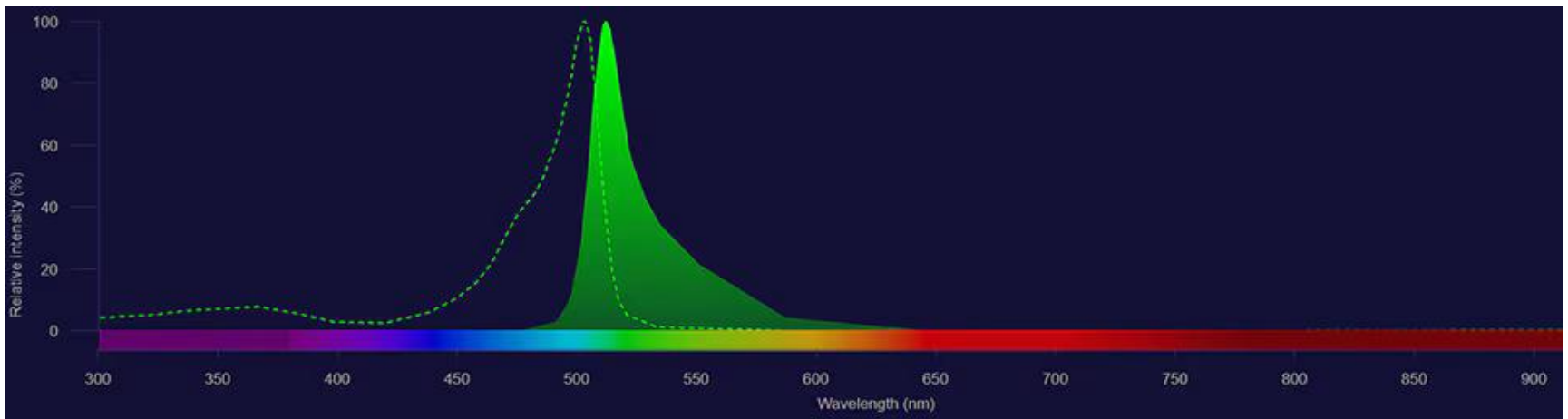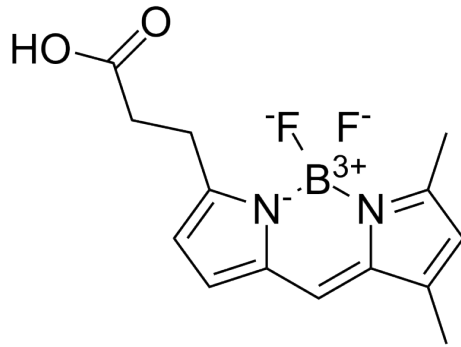

Announce

Concepts

Procedure



Hazards

Tips



Clarification

End

Intro to BODIPY-FL

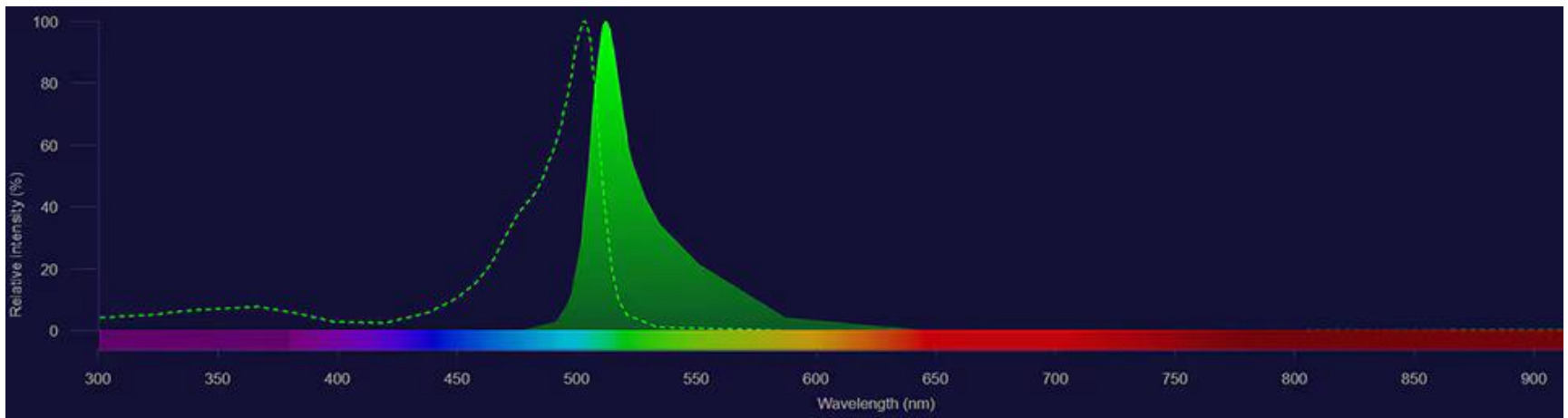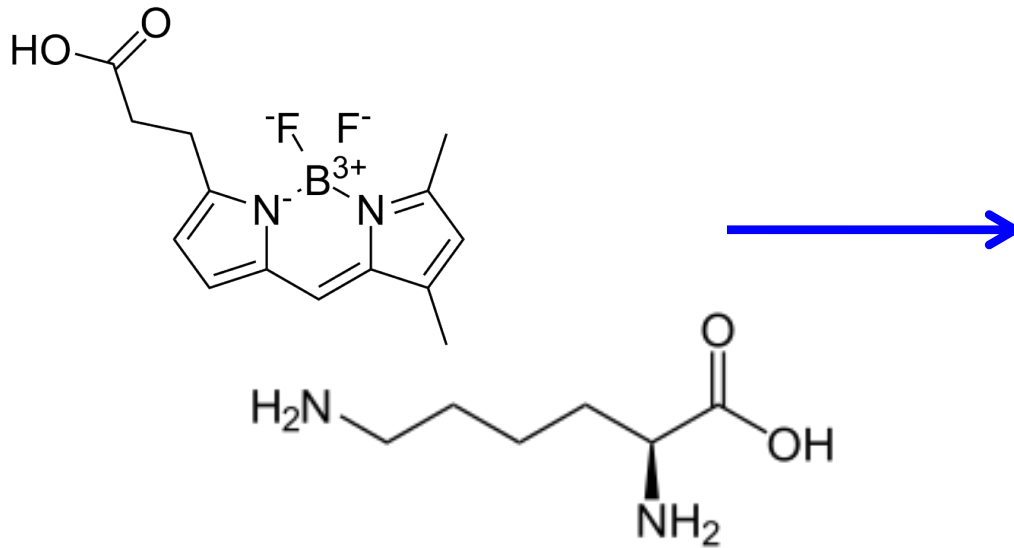

Intro to BODIPY-FL

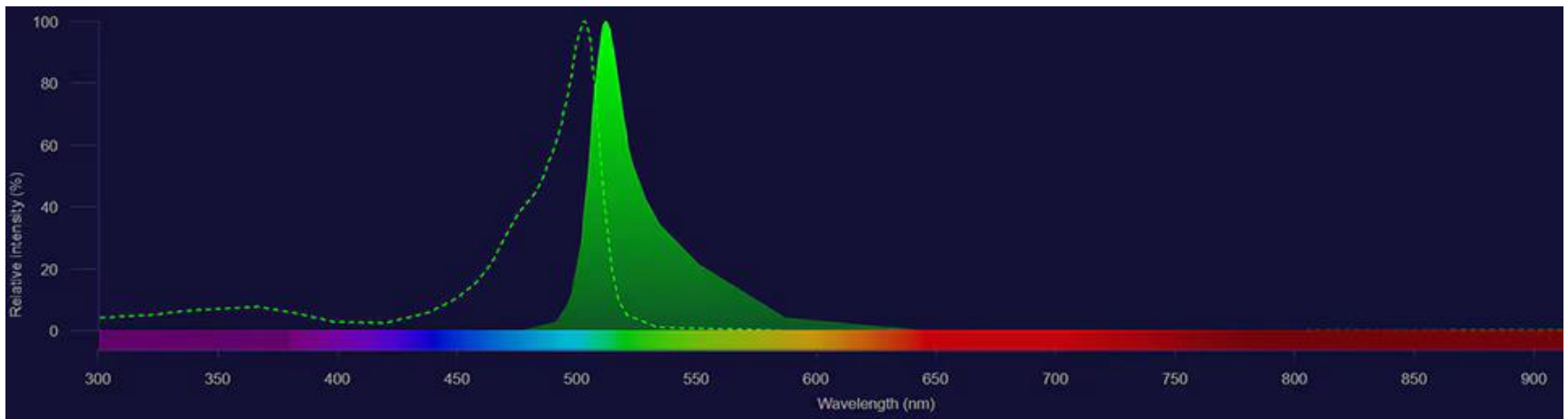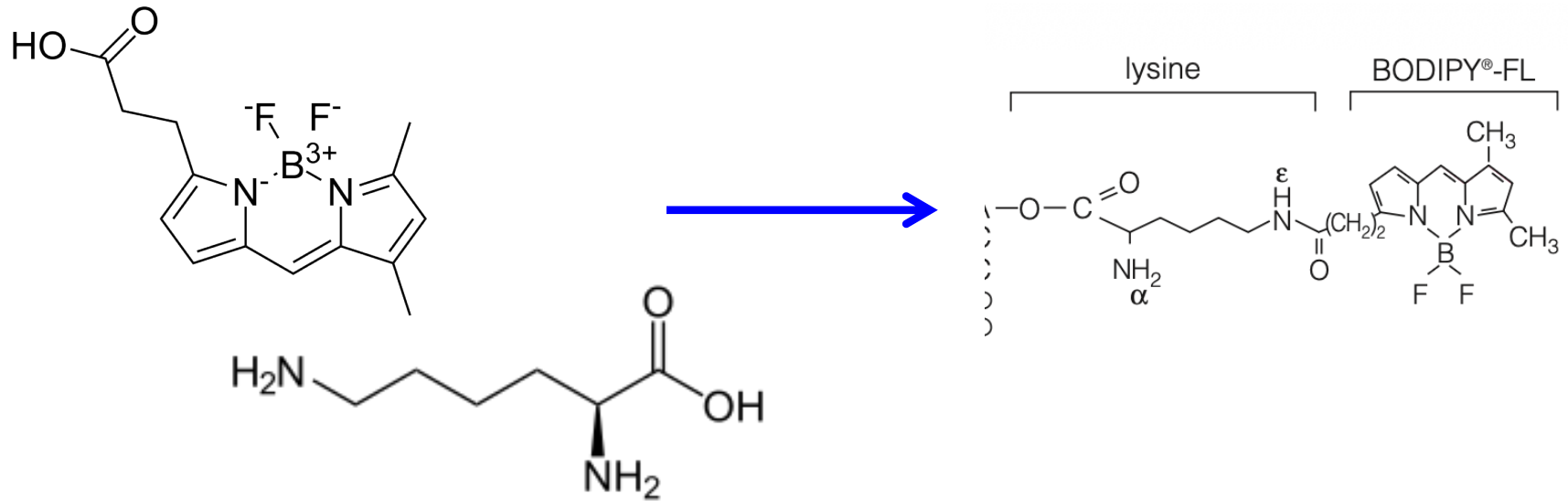

Nothing is forever

- Fluorophores can only fluoresce a limited number of times
- If they are under room light for too long, you won't be able to see them on the gel.
- When samples are in the ice bucket, cover with lid.
- When covering with the bucket lid is not practical, wrap in aluminum foil.

Properties of RNases

- RNA is easily degraded by ribonucleases, RNases
- These enzymes are on your clothes, skin, saliva
- Very important to work with gloves
- You can inactivate RNases by sterilizing solutions at high temperatures in the presence of DEPC, or by using chemical modifications

Properties of RNases

- RNA is easily degraded by ribonucleases, RNases
- These enzymes are on your clothes, skin, saliva
- Very important to work with gloves
- You can inactivate RNases by sterilizing solutions at high temperatures in the presence of DEPC, or by using chemical modifications

We want to make RNA transcripts, be careful of RNase contamination!

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
Chapter 8AB: Procedure

- Use gloves at **all times** to prevent RNase contamination
- We will provide sterile/RNase-free pipet tips, centrifuge tubes, etc.
Please keep them that way!!!
- You will be pipetting **very small volumes!**
 - We should have some P-2 pipets available (**0.2 – 2 μ L**)

Chapter 8AB: Procedure

- Tips for working with small volumes:
 - Spin down samples in tabletop centrifuge
 - Pipette samples as little as possible
 - Pipette larger volumes into smaller volumes
 - Check pipette tips

Chapter 8AB: Reagent List (example)– make your own! It will be different from this

<u>Reagent Name</u>	<u>Cap mark</u>
Nuclease-free water	
TNT reaction buffer	
Amino Acid mix –Met	
RNasin RNase Inhib.	
TNT SP6 RNA polymerase	
Fluortect GreenLys	
TNT Wheat Germ Extract	
SP6-Luciferase (0.5 µg/µL)	
pGEM3-Rel (0.5 µg/µL)	
pGEM4-Rel (0.5 µg/µL)	

Chapter 8AB: Procedure

A. Transcription/Translation:

- Work together with a second group for the first part
- Thaw reagents
- Set up cocktail for 4 reactions according to protocol
- Prepare mastermix in tube pre-aliquoted with your fluorescent lysine
- Remove 5 μL x 2 for zero-time point control reactions
- Aliquot cocktail, 19 μL / reaction and add 1 μL DNA to each
 - pGEM3-Rel
 - pGEM4-Rel
 - Positive control – DNA template encoding luciferase (SP6-luciferase)
 - Negative control – No DNA (nuclease-free water)

Chapter 8AB: Procedure

A. Transcription/Translation:

- Incubate all samples at **30 °C** for **90 min**
- Stop reactions on ice
- Split reactions – 10 μ L/group
- **Each group works separately from now on**
- Prep samples for SDS-Page gel
- Store samples at -20 °C until next lab

Chapter 8AB: Procedure

- You will be supplied with the positive control plasmid “SP6-Luciferase” containing the *luciferase* gene
 - This plasmid has the optimal coding sequence for the SP6 promoter
 - Lots of RNA > lots of protein > a lot of fluorescence
- You may choose to use your own A & B plasmids (check concentration and determine how much to use)
- We can also supply you with A & B plasmids

Chapter 8AB requirements

- Please make sure to submit your Pre-Lab to GradeScope *before* your lab section
- Include in your procedures:
 - a flowchart for the transcription/translation protocol
 - table for the code for reagent tubes
- Please submit your in-lab data collection assignment to Gradescope *by the end of the day after* your lab section

Chapter 8AB

Before the lab period, you should have:

- ✓ Completed your Pre-lab Write-up and submit on Gradescope
 - ✓ Title, purpose and procedures
 - ✓ Remember to include:

At the end of lab, you should have:

- ✓ Performed transcription/translation on 4 samples
- ✓ Saved SDS-PAGE samples for next week (the bulk of the radioactivity!).
Be sure to label well.

Questions?

**In-class activity
&
Discussion Quiz**